# Aseptic sampling valve

## Type VPA

### CHARACTERISTICS

The AR sampling valve range has been designed for sterile sampling applications in the Chemical and Pharmaceutical industries. Its compact design allows easy installation on small vessels or process lines.

### FULLY AUTOCLAVABLE,

Valve body and head are made out of certified 316L bar stock.

Damage to the seating plug and unexpeted material in the outlet are indications of overpressure. A leak detector hole is placed immediately behind the outlet. On request the valve can be equipped with a CIP system.

### SURFACE FINISH

External mechanical "mirror" polish  
Internal mechanical polish  $Ra < 0.4 \mu m$

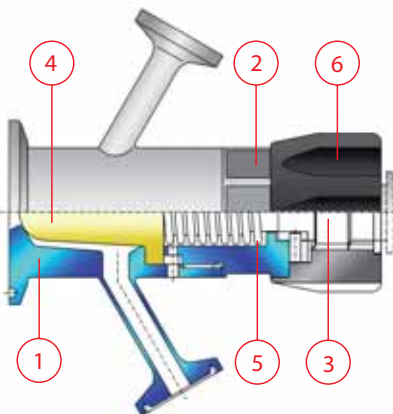
### VALVE BODY connections available in:

- weld ends
- clamp ends

Other types available upon request.

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

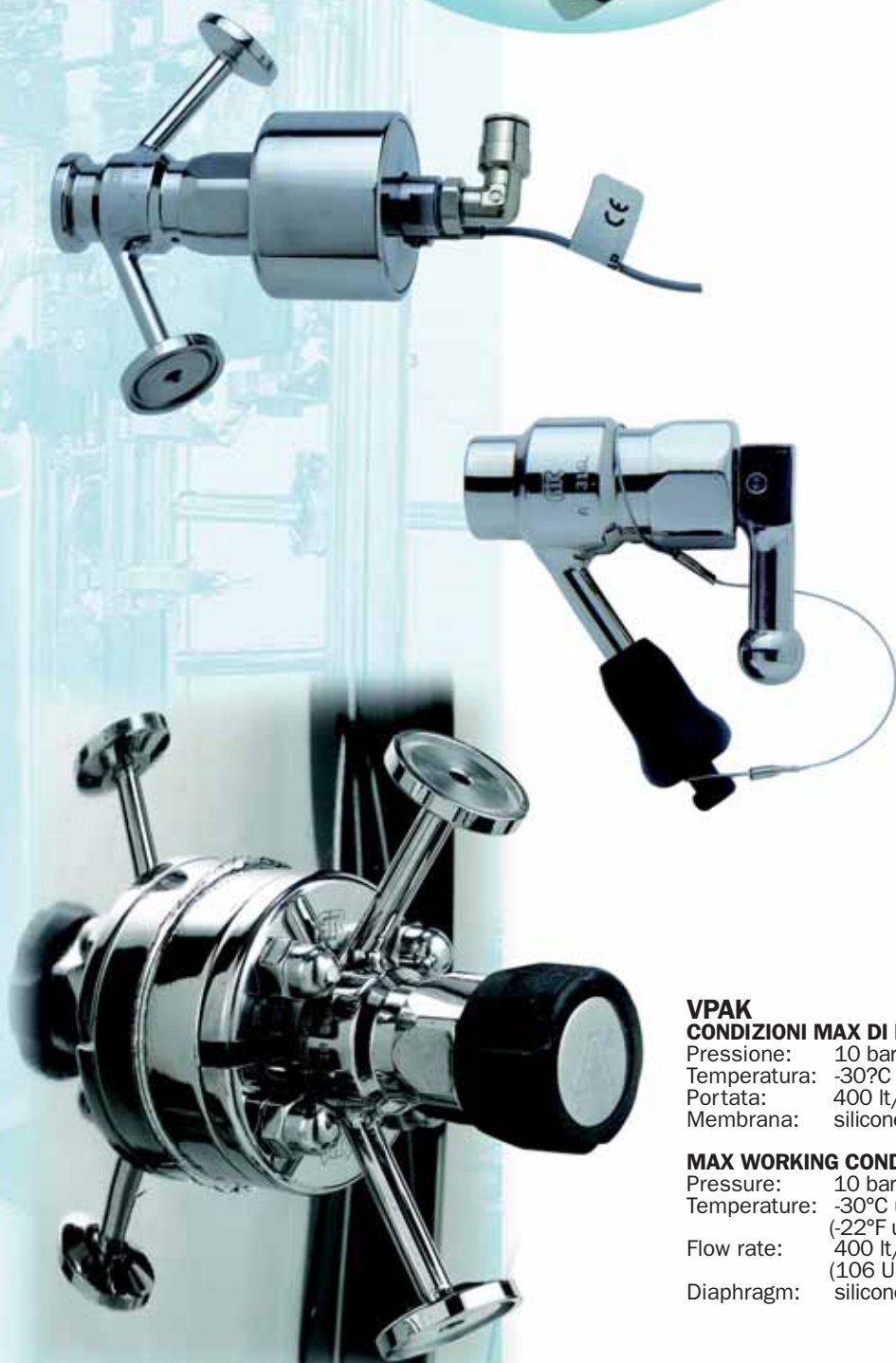
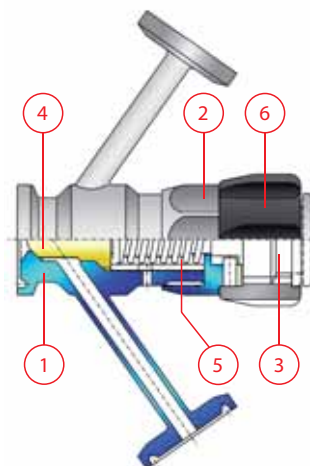
## STERILE SAMPLING

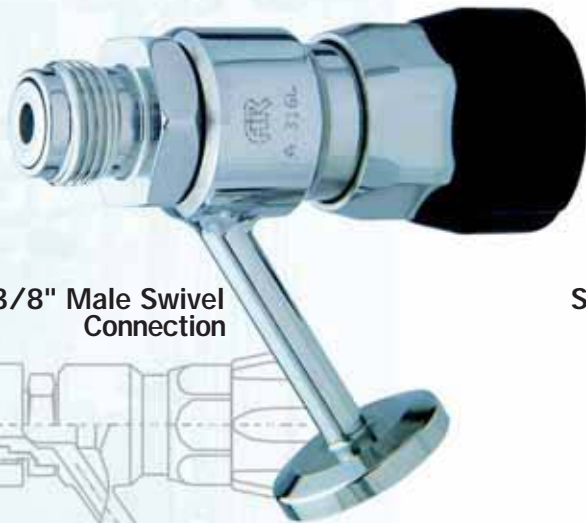Developed for taking samples in a sterile way, the special configuration assures effective cleaning (CIP) and sterilization (with saturated steam) of product contact surfaces of valve and seating plug with the valve in closed position. Once installed, the sealing plug shuts off flush with the inlet port connection with a zero dead leg.

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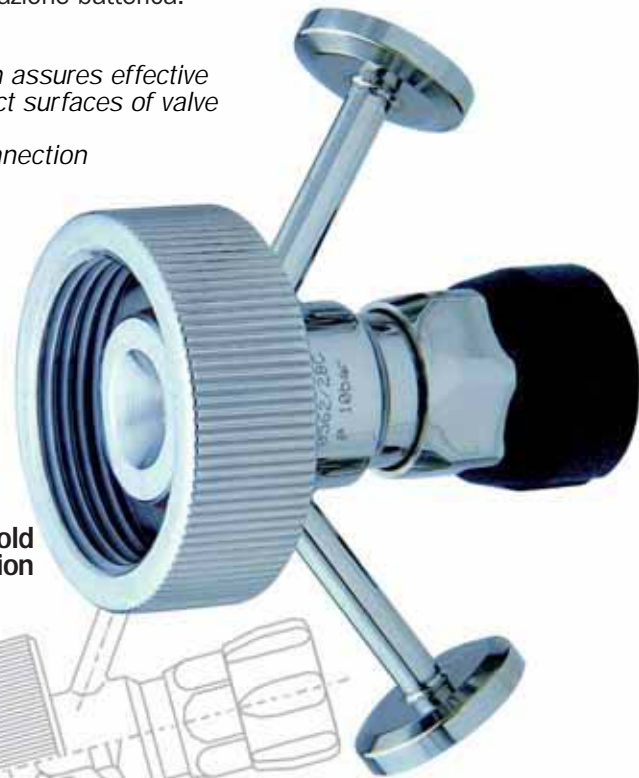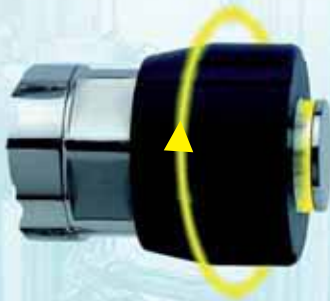
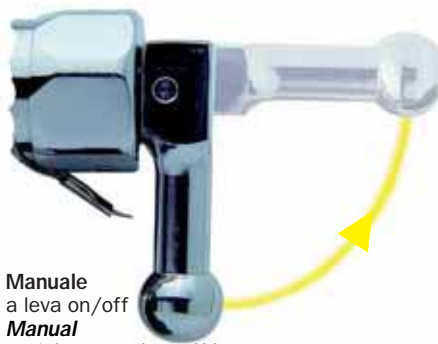
**Manuale**  
a leva on/off  
**Manual**  
stainless steel on-off lever

**Pneumatico** inox on/off  
**Air operated** stainless steel air-spring actuators

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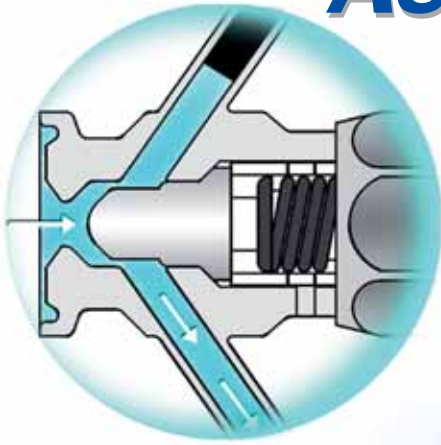
## DOCUMENTATION

All valves are fully traceable for validation processes. Each valve is accompanied by material test certificates EN 10 204 3.1B 3A standards compliance certificates, FDA regulations compliance certificates, and tapes detailing the valve interior surfaces Ra.

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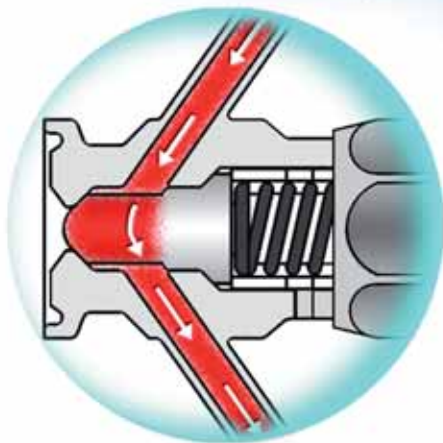
### IMPORTANT!

For reliable performance there must be a safety margin between the spring load of the seating plug and the maximum pressure in the product line.

Higher pressures will force open the valve.

Please indicate the maximum pressure of your applications.

## SIP phase

■ Vapore  
Steam

**fig. 2**  
Valvola chiusa (fase di CIP/steriliz.)  
Valve closed (CIP/steril. phase)

## Come ordinare How to order

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**1**   **2**   **3**   **4**   **5**   **6**  
**VPAK / M / ETC / UTC / VTCX**

Tipo valvola  
Valve type  
VPA  
VPAK

Attuatore  
Actuator  
M Manuale  
Manual knob  
ML Manuale leva  
Manual lever  
PN Pneumatico  
Pneumatic  
PS Pneum + Switch

Attacco corpo  
Valve connectors  
ETC clamp BS  
ESD a saldare  
weld  
EGM filettato  
threaded  
EFL flangiata  
flanged  
EIN ingold  
connection

Attacco sample  
Sampling connectors  
UTC clamp BS  
USD a saldare  
weld  
UPG a portagomma  
rubber

Attacco vapore  
Steam connectors  
VTC clamp BS  
VSD a saldare  
weld  
VPG a portagomma  
rubber

Membrana  
Diaphragm  
X silicone  
G PTFE