# Aseptic Sampling Valves

Sanitary  
flow  
equipment

Tipo **VPK**

# Presacampione aseptica

## CARATTERISTICHE

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## STERILE SAMPLING

Developed for taking samples in a sterile way, the special configuration assures effective cleaning (CIP) and sterilization (with saturated steam) of product contact surfaces of valve and seating plug with the valve in closed position. Once installed, the sealing plug shuts off flush with the inlet port connection with a zero dead leg.

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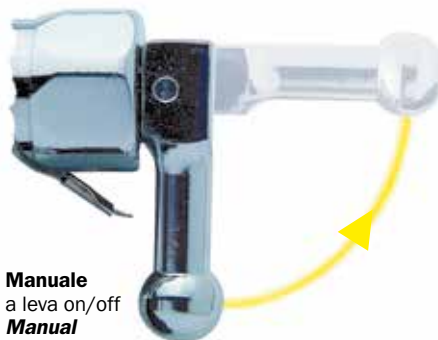
**Manuale**  
a leva on/off  
**Manual**  
stainless steel on-off lever

**Pneumatico** inox on/off  
**Air operated** stainless steel air-spring actuators

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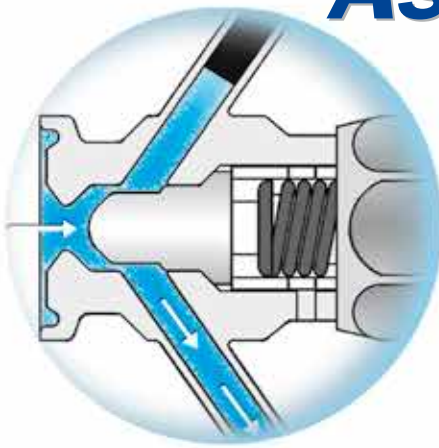
## DOCUMENTATION

All valves are fully traceable for validation processes. Each valve is accompanied by material test certificates EN 10204 3.1, FDA regulations compliance certificates, and surface roughness conformity certificates.

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
## IMPORTANT!

For reliable performance there must be a safety margin between the spring load of the seating plug and the maximum pressure in the product line. Higher pressures will force open the valve. Please indicate the maximum pressure of your applications.

## Sampling phase

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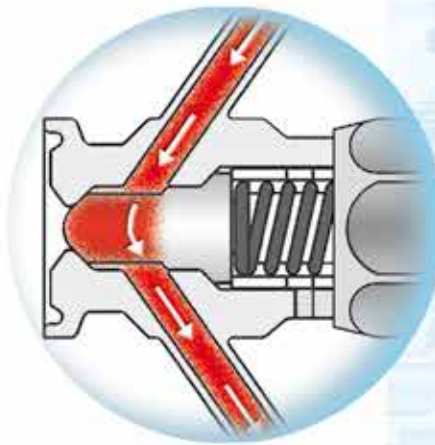
**fig. 1**  
Valvola aperta (fase di prelievo)  
Valve open (sampling phase)

## CIP/SIP phase

■ Vapore  
Steam

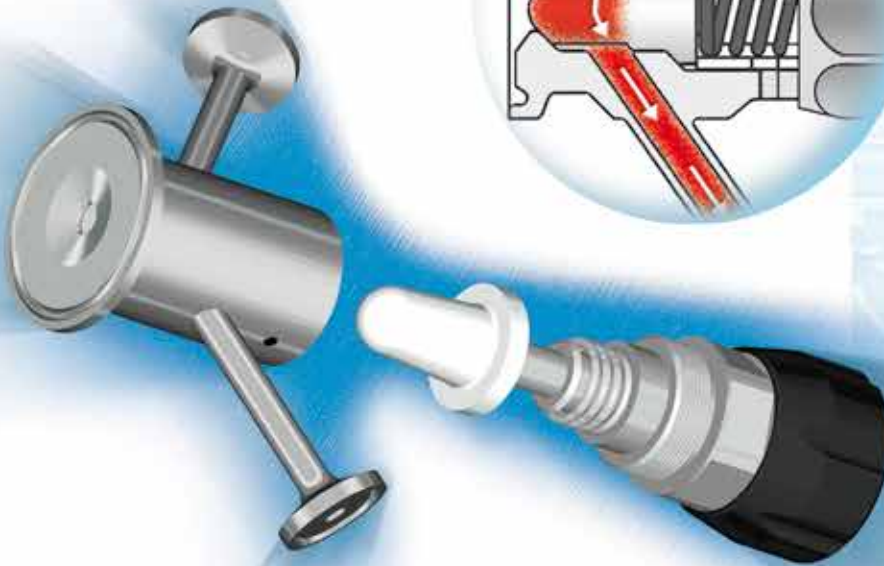**fig. 2**  
Valvola chiusa (fase di CIP/steriliz.)  
Valve closed (CIP/steril. phase)

## VPA - VPK ORDERING INFORMATION

To specify the part completely, start with the product description and select the additional options as shown below:

Valve	Option	Inlet hole	Inlet	Sampling	CIP/SIP	Diaphragm	Actuator	Sensor	finish	material
VPK	-	05	AG	PB	AG	T	MC		-11	A

Valve	Inlet Hole	Inlet Connection	Sample - CIP/SIP	Diaphragm	Surface Finish VPK
VPK	05 5 mm	1/2" AG •	clamp 1/2" AG • •	T TFM PTFE	11 ≤ 0,5 µm
VPA	07/10/13 7 / 9,4 / 12,7 mm	3/4" AH • •	clamp 3/4" AH • •	X Silicone	51 ≤ 0,5 µm + Electropolish
VPZ	05/07/10/13	1" AJ • • •	to be weld Ø6,25x1,24mm RB • •		Surface Finish VPA
on request (*) 7mm Silicone only		1 1/2" AM • • •	to be weld Ø12,7x1,65mm BG • •		71 ≤ 0,5 µm
		2" AN • • •	hose barb Ø6,25mm PB • •		41 ≤ 0,4 µm + Electropolish
		butt weld Ø19mm QC • •	hose barb Ø12,7mm PF • •		on request
		Ø30mm QF • • •	quick connect ZG • • •		
		1/4" LC • • •	on request		
		BSP screw 3/8" LE • • •			
		1/2" LG • • •			
		3/4" LH • • •			
		Ingold Ø25 ZA • • •			
		flanged S... • • •			
		on request			

Option	Actuator	Sensor	Materials
- Standard Safe Supply	MC Manual wheel	11 Inductive PNP 12/24VdC	A AISI 316L-1.4404
A ATEX execution	ML Manual lever on/off	on request	B 1.4435
	MR Lever auto return		C C22 Hastelloy
	PN Pneumatic s.e./n.c.		D C276 Hastelloy
	PL Pneumatic + lever		on request

# Aseptic sampling valve

## Type VPA

### CHARACTERISTICS

The AR aseptic sampling valve range has been designed for sterile sampling applications in the Chemical and Pharmaceutical industries. Its compact design allows easy installation on small vessels or process lines.

### FULLY AUTOCLAVABLE,

Valve body and head are made out of certified 316L bar stock.

The aseptic design ensures that, during sampling, the fluid enters into contact only with the valve body and the diaphragm, excluding the mechanical parts of the actuator from the process.

Upon request they are equipped with a specular connection to the sample that allows washing by CIP or sterilization by SIP. Available with manual or pneumatic actuator that are perfectly interchangeable with each other.

The valve body is equipped with a leak hole that indicates the state of the membrane.

### SURFACE FINISH

External mechanical "mirror" polish

Internal mechanical polish  $Ra \leq 0.5 \mu m$

### VALVE BODY connections available in:

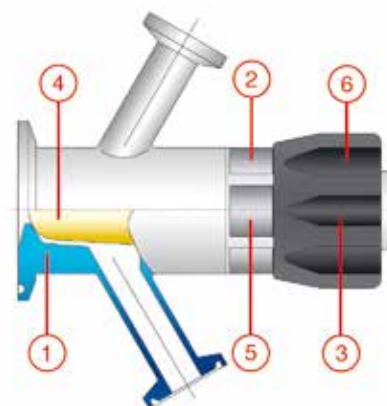
- weld ends

- clamp BS 4825 / ASME BPE

Other types available upon request.

### MATERIALS

- 1 AISI 316L valve body
- 2 AISI 316L valve head
- 3 AISI 316L stem
- 4 Silicone or PTFE membrane
- 5 AISI 304 spring
- 6 adjustable autoclavable PTFE knob

### VPA

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Pressione: 10 bar (g)  
Temperatura: +150 °C  
Portata: 1.200 lt/h a 2 bar  
Membrana: silicone - PTFE

#### MAX WORKING CONDITIONS

Pressure: 10 bar (145 Psi)  
Temperature: +150 °C  
+302 °F  
Flow rate: 1.200 lt/h at 2 bar  
(317 US gal/hr at 29 Psi)  
Diaphragm: silicone - PTFE

ASEPTIC SAMPLING VALVES

SANITARY SAMPLING VALVES

SANITARY SAMPLING BOTTLE

TANK BOTTOM ASEPTIC VALVE

SPRING CHECK VALVES

HIGH PURITY BALL VALVES

BUTTERFLY VALVES

DTS HEAT EXCHANGERS

SIGHT GLASS-FLOW INDICATOR

CLAMP FITTINGS

TANK CONNECTIONS

SILICONE HOSE & FITTINGS

MAGNETIC MIXER

ITT DIAPHRAGM VALVES

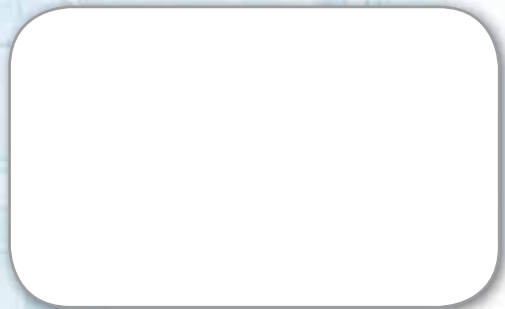
WASHING DEVICES

Sanitary flow equipment

AERRE INOX s.r.l.  
Via Gerola, 4  
I-26010 Fiesco (Cremona)  
Tel. +39 0374 370 828 Fax +39 0374 370 833  
www.aerreinox.it

Get the information you need and more at [info@erreinox.it](mailto:info@erreinox.it)